Tris-NaCl buffer

REF ARO04K

Bf 3 flacons x 50 mL / 3 vials x 50 mL

Tris-NaCl buffer
FOR RESEARCH USE ONLY.
NOT FOR USE IN DIAGNOSTIC PROCEDURES

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English, revision: 09-2022

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INTENDED USE:

Dilution buffer for coagulation assays, or for any application where Tris Buffer Saline (TBS) is needed. Ready to use.
This kit is for research use only and should not be used for patient diagnosis or treatment.

REAGENTS:

Bf **Tris-NaCl buffer**, pH : 7.50, Ionic Strength: I=0.15, liquid form. Contains Tris at approximately 0.05M (Tris (hydroxymethyl) aminomethane), Sodium Chloride at approximately 0.15 M, preservatives and stabilizers.

The product is classified as non-hazardous and is not subject to labeling according to EC Regulation No. 1272/2008 [CLP].

WARNING AND PRECAUTIONS:

- Waste should be disposed of in accordance with applicable local regulations.
- Use only the reagents from the same batch of kits.
- This device of *in vitro* use is intended for professional use in the laboratory.

REAGENT PREPARATION:

Bf Reagent is ready to use, allow to stabilize for 30 minutes at room temperature (18-25°C), homogenize before use.

STORAGE AND STABILITY:

Unopened reagents should be stored at 2-8°C in their original packaging. Under these conditions, they can be used until the expiry date printed on the kit.

Bf Reagent stability after opening, free from any contamination or evaporation, and stored closed, is of:

- 4 weeks** at 2-8°C.

LIMITATIONS:

- Any reagent presenting no limpid appearance or showing signs of contamination must be rejected.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

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